




Perfusion-Set for "60" Petri dishes

500-800 260

Perfusion-Set for "60" Petri dishes

consisting of:

cover with glass insert, # 000 861; perfusion ring, # 501 466;
canal tube inlet, # 501 425; canal tube outlet, # 501 426;
grub screws # 900 094; screwdriver # 900 405

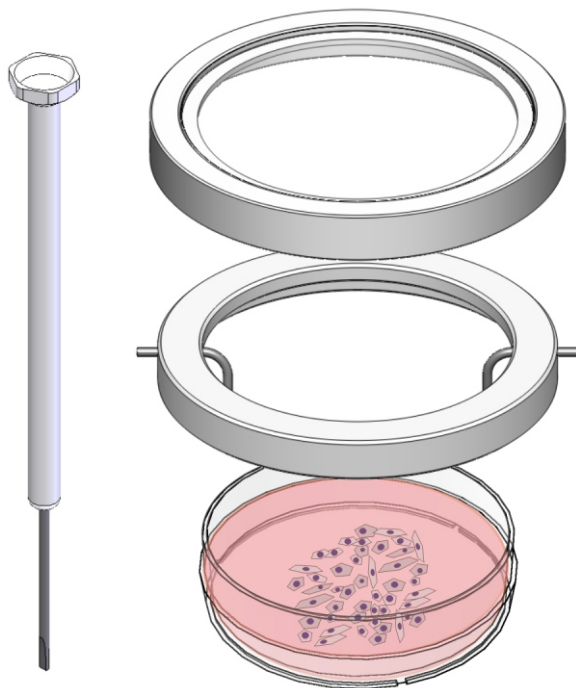

cover with glass insert

- stainless steel V4A
- sterilizable
- glass insert
- cut-outs for the canal tubes

perfusion ring


- stainless steel
- sterilizable
- lateral openings for the canal tubes
- Screws for the fixation of the canal tubes

Petri dishes "60"

- different manufacturer with outer diameter 50,5 - 56,0 mm

Scheme of an "open perfusion"

assembly and procedure of the perfusion

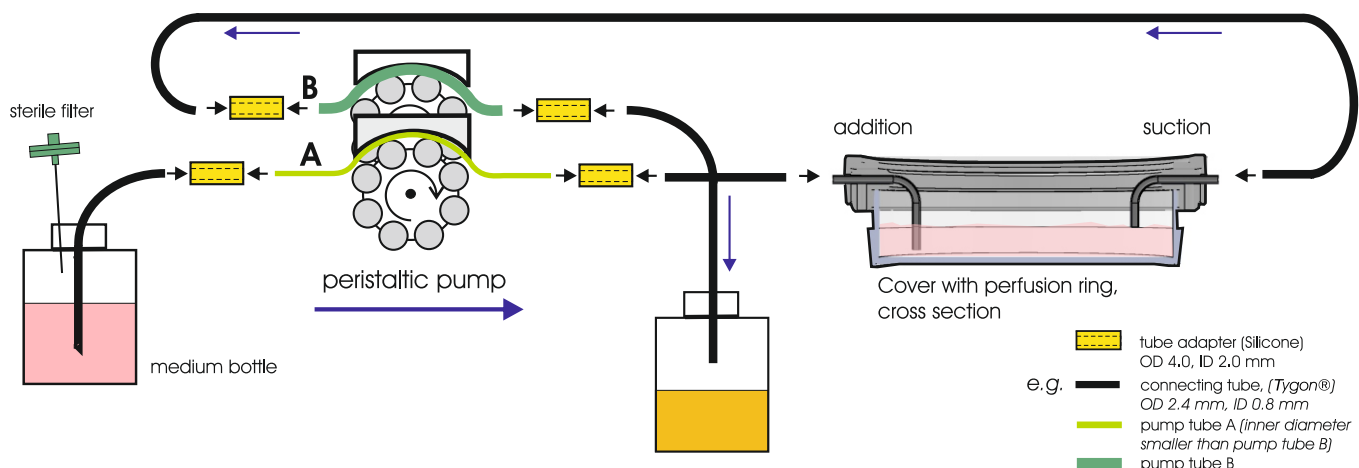


This Perfusion-Set is compatible to a large variety "60" Petri dishes.

**Please do not hesitate to contact us regarding a
cover with perfusion adapter for other Petri dishes!**


Perfusion-Set for "60" Petri dishes

500-800 260

cover with glass insert

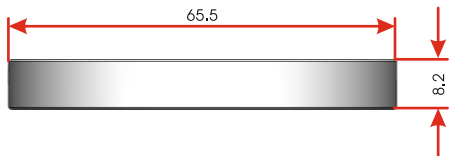


perfusion ring with canal tubes

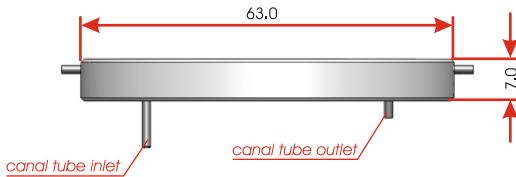


Specifications

perfusion ring with cover - bottom view

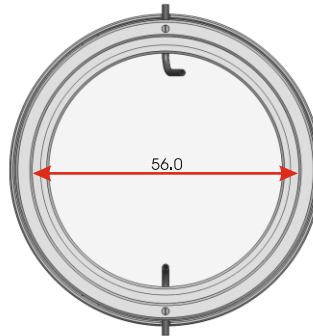

dimensions in mm

- For "60" Petri dishes with an outer diameter:
Dmax = 56.0 mm
Dmin = 50.5 mm

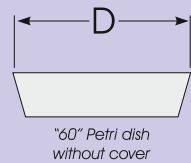



Assembly

- Set the sterilized perfusion ring onto the Petri dish "60"
- Set the cover with glass insert onto the perfusion ring
- Plug the Tygon[®] tubes onto the canal tubes



Application example

- Petri dish "60" with perfusion ring and cover inserted into the Heating Insert P 2000

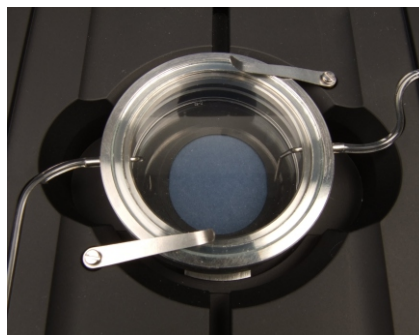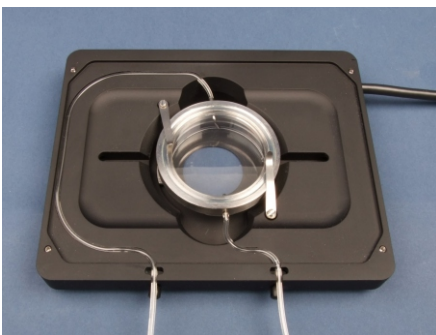

This Perfusion-Set is compatible to a large variety "60" Petri dishes.

Please do not hesitate to contact us regarding a cover with perfusion adapter for other Petri dishes!