


## Perfusion-Set for "60" Petri dishes

# 500-800 260

### Perfusion-Set for "60" Petri dishes

consisting of:

cover with glass insert, # 000 861; perfusion ring, # 501 466;  
canal tube inlet, # 501 425; canal tube outlet, # 501 426;  
grub screws # 900 820; screwdriver # 900 405

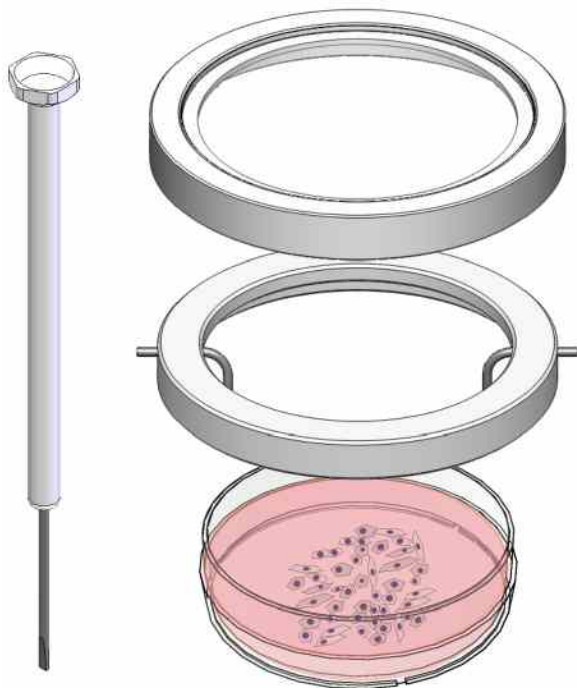

cover with glass insert

perfusion ring

Petri dishes "60"


- stainless steel
- sterilizable
- glass insert
- cut-outs for the canal tubes

- stainless steel
- sterilizable
- lateral openings for the canal tubes
- Screws for the fixation of the canal tubes

- different manufacturer with outer diameter 50.5 - 56.0 mm

### Scheme of an "open perfusion"

assembly and procedure of the perfusion

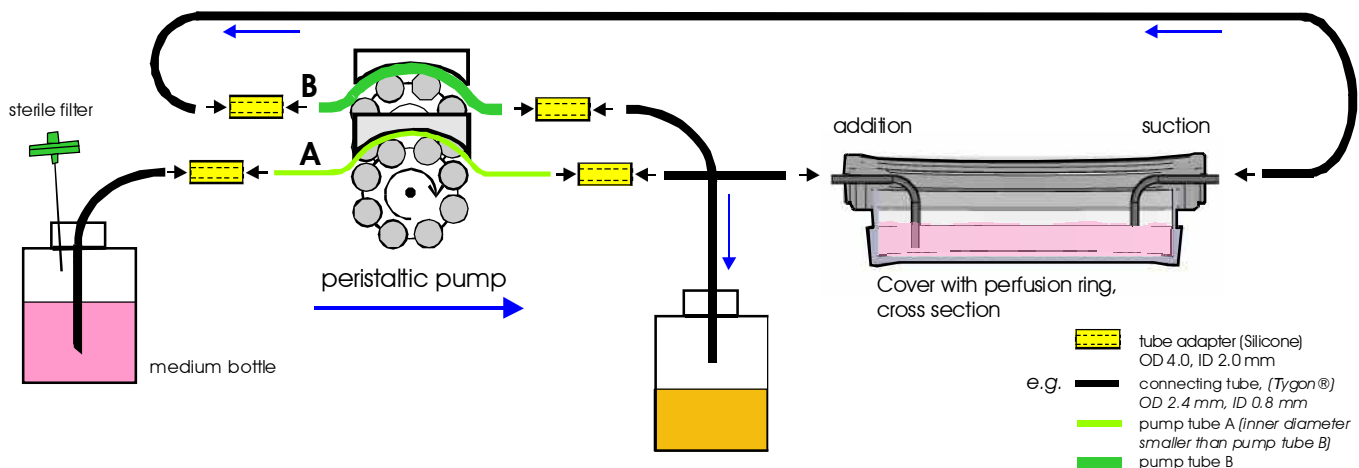

This Perfusion-Set is compatible to a large variety "60" Petri dishes.


**Please do not hesitate to contact us regarding a cover with perfusion adapter for other Petri dishes!**

page 1/2

Version 1.1, 10-2017  
[www.pecon.biz](http://www.pecon.biz)

## Perfusion-Set for "60" Petri dishes

# 500-800 260

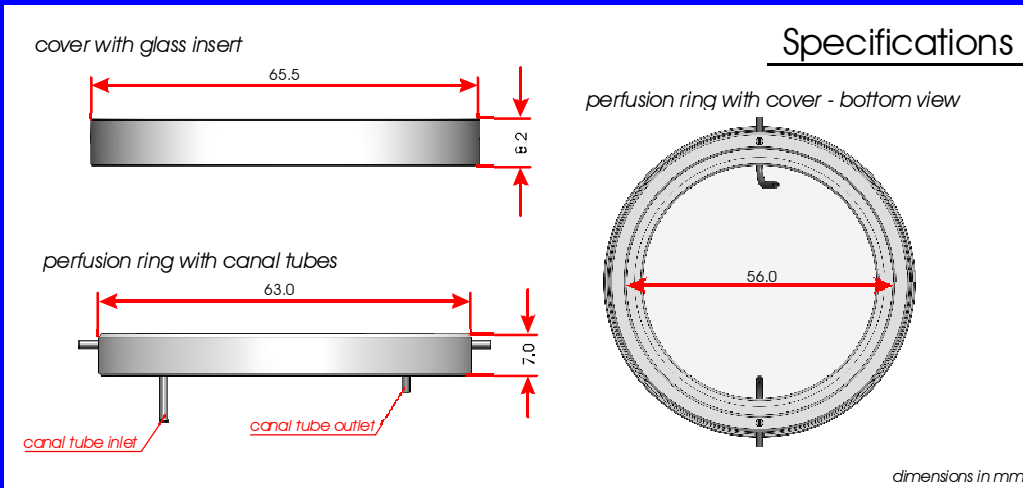

### Specifications


- For "60" Petri dishes with an outer diameter:  
Dmax = 56.0 mm  
Dmin = 50.5 mm

### Assembly



- Set the sterilized perfusion ring onto the Petri dish "60"
- Set the cover with glass insert onto the perfusion ring
- Plug the Tygon<sup>®</sup> tubes onto the canal tubes

### Application example

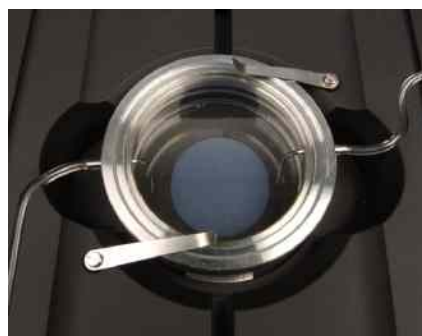

- Petri dish "60" with perfusion ring and cover inserted into the Heating Insert P 2000

This Perfusion-Set is compatible to a large variety "60" Petri dishes.  
**Please do not hesitate to contact us regarding a cover with perfusion adapter for other Petri dishes!**